

Do not fill 96 well plates over “2/3” of maximum volume, you will lose significant amount of your sample by overfilling.

Non-Skirted Plates need adaptor in drawer below

Maximum weight per plate and adaptor is 115 grams.

Load plate samples symmetrically see diagram below.

Concentrator setting is V-AQ.

Brake ON

Heat according to sample size, this will vary.

Time according to sample size, this will vary.

Manually vent chamber during run by pressing “mode vent” at middle of run time and at the end of run time to clear moisture from tubing.

Clean any spills immediately to avoid staining chamber.

Leave lid open when finished.

Turn off.